

DfCell Mycoplasma Off Spray

Product description

DfCell Mycoplasma Off Spray effectively inhibits the growth and contamination of fungi (including spores), bacteria (including spores), mycoplasma, and viruses (including HIV and Hepatitis B). Its active ingredient is a quaternary ammonium derivative. It is free from mercury, formaldehyde, phenol, alcohol, and other irritants, non-toxic, and biodegradable. It does not cause any damage to general work surfaces and equipment.

This product is a ready-to-use solution, to be used directly.

Specifications

Catalog Number	C230101E/C230101S/C230101M
Specifications	500 mL /2×500 mL/10×500 mL

Storage

Store and transport at room temperature. Shelf life is 18 months.

Notes

1. Please read the entire instruction manual carefully before using this reagent.
2. Standardized operations, including the preparation of the reaction system, sample handling, and sample addition, should be followed throughout the experiment.
3. For your safety and health, wear laboratory coats and disposable gloves during operation.
4. For research use only.

Instructions

1. Spray the CO₂ incubator once every two weeks. There is no need to empty the incubator before spraying.
2. Spray the laminar flow hood before each use and allow it to air dry naturally.