

ATP Luminescent Cell Viability Assay Kit

Product Information

Product name	Cat#	Specification
ATP Luminescent Cell Viability Assay Kit	40210ES10	10 mL
	40210ES60	100 mL
	40210ES80	10×100 mL

Product Description

ATP is the most important energy molecule in cells, which can be used to measure the cell metabolism level and has a good linear relationship with the number of living cells. Therefore, the number of living cells can be reflected by ATP content. This kit uses the chemiluminescence signal to measure ATP content in cells through the luciferin luminescence reaction catalyzed by ATP-dependent luciferase, so as to detect cell viability or quantitative detection of the number of living cells, with high sensitivity and wide linear range. This kit is compatible with a small number of samples and a large number of high-throughput screening tests.

The luminescent cell viability detection reagent provided by this kit has a wide linear range, high sensitivity and good stability. In the 96-well plates, there was a good linear relationship between 100 and 100,000 cells, but the upper limit for the number of cells was different. In addition, the operation is simple, the detection reagent provided in the kit is ready-to-use, stable reading, fast detection, only about 10 minutes to complete the detection, there is no need to wash cells, and there is no need to replace or remove the culture medium. Compared with other common cell viability determination methods, such as Calcein-AM, CCK-8, etc., the cell viability detection by luminescence method is simpler and faster.

Shipping and Storage

The components are shipped with ice pack and can be stored at -20°C for 1 year.

Cautions

1. The reagent contains luciferase, and repeated freezing and thawing will affect its activity. It is recommended to store at -20°C away from light.

2. Reagents and cell samples should be balanced to room temperature before use to avoid the influence of enzyme catalytic effect.

3. High drug content may interfere with luciferase reaction, thus affecting chemiluminescence signal. It is recommended to set up cell culture medium containing drugs to illuminate the hole to eliminate solvent interference.

4. For your safety and health, please wear lab coat and disposable gloves.

5. For research use only!

Instructions

1. Cell culture

A 96-well plate suitable for chemiluminescence detection was used, and 100 μ L cells were inoculated in each well (the initial inoculated cell density was determined according to the culture time, and the number of cells in each well should not exceed 100,000). Meanwhile, the Wells without cells were set as negative control. Cells were cultured at 37°C with 5% CO₂. Cell concentration gradients can also be set to obtain optimal experimental results. Treat the cells at appropriate times as needed.

2. (Optional) ATP standard curve production

Dilute the prepared ATP standard solution to an appropriate concentration gradient with PBS, add 100 μ L standard solution to each well of 96-well plate.

3. Cell viability detection

YEASEN

1. Melt and thaw the existing luminescence detection reagent, and balance it to room temperature;

2. Take out the cell culture plate and balance it at room temperature for 10 mins;

3. Add 100 µL detection reagent to each well of 96-well plate (due to edge effect of well, it may lead to unstable luminous signal, so

it is not recommended to lay plate at the edge);

4. Oscillate at room temperature for 2 mins to promote cell lysis;

5. Place at room temperature for 10 mins to stabilize the luminescent signal;

6. Chemiluminescence detection was carried out with multi-functional enzyme reader. Set the corresponding parameters according to the requirements of the instrument, the detection time of each hole is generally 0.25-1 s, the specific need to be adjusted according to the sensitivity of the instrument;

7. Calculate the relative activity of cells according to the chemiluminescence reading, or calculate the ATP content according to the ATP standard curve to obtain the relative activity of cells.

[Notes] : The detection effect varies with the type of cells. For some cells with particularly high ATP content, chemiluminescence readings may continue to rise but lose linear relationship when the number of cells reaches 100,000 or more.