

Spray (Ready-to-use)

Product Information

Product Name	Cat#	Specification
	40605ES02	500 mL
Spray (Ready-to-use)	40605ES03	2×500 mL
	40605ES08	10×500 mL

Product Description

Spray (Ready-to-use) stops the growth and contamination of fungi (and spores), bacteria (and spores), mycoplasmas and viruses (including HIV and Hepatitis B). Its active ingredients are quaternary ammonium derivatives, free of irritating substances such as mercury, formaldehyde, phenol, or alcohol, non-toxic and biodegradable. There is no damage to the surface of the general workbench and equipment.

This product is a working solution of 1×, which can be used directly.

Instructions

- 1) Spray the CO₂ incubator once every two weeks, no need to empty the incubator before spraying.
- 2) Spray the ultra-clean workbench once before each use and let it dry naturally.

Shipping and Storage

Shipping and store at room temperature, the shelf life is 18 months.

Cautions

- 1) Please read this manual carefully before using this reagent.
- 2) The entire experiment should be standardized, including the preparation of the reaction system, sample processing and sample addition.
- 3) For your safety and health, please wear a lab coat and disposable gloves for operation.
- 4) For research use only!